



Note: Read the corresponding technical documentation for handling and safety reasons.

Science with Passion



AZURA® SmartClick StartUp Kit Installation instructions

The AZURA SmartClick StartUp Kit includes two SmartClick Edges (one for the back of the device and one for the side) and a SmartClick Rail. AZURA SmartClick works only with AZURA L devices. You can attach SmartClick Edges to both sides of an AZURA L device. Up to two SmartClick Rails can be mounted on the SmartClick Edges. The number of modules on a SmartClick Rail depends on their size, especially when using a second rail.

These instructions explain how to install the SmartClick Edges for the SmartClick Rail(s).

Tools required
(included in accessories)

- Screwdriver X0257 (Torx T10)
- Screwdriver X0010-1 (hexagon socket) for older devices

Procedure

Process

1. On the back of the device, remove the two screws ② that secure the side panel ① to the housing (shown in the illustration on the left side of the device).

Figure



Fig. 1: Removing the two screws on the back of the device

2. On the same side of the device, attach the SmartClick Edge ③ to the side of the device using the screwdriver.

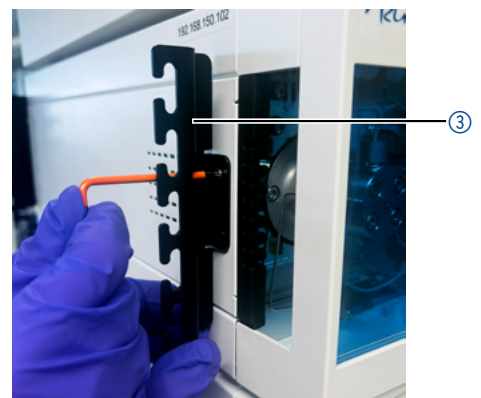


Fig. 2: Attaching the SmartClick Edge to the side of the device



Note: First, attach the SmartClick Edge to the side. If you attach the SmartClick Edge to the back of the appliance first, the holes on the side will shift so that you will no longer be able to screw the screws in.

Process

3. Attach the second SmartClick Edge ④ to the back of the device using the screwdriver.



Note: Use the screws from the accessory pack to install the SmartClick Edges. The screws that are originally installed are not long enough.

Figure

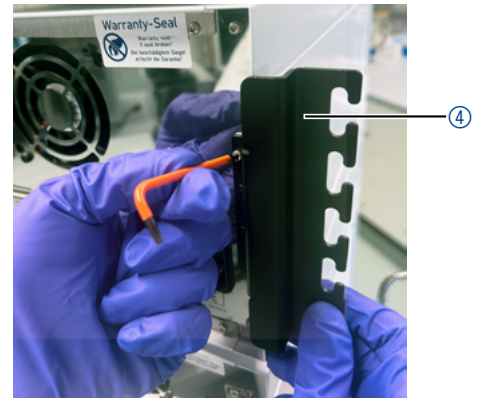


Fig.3: Attaching the SmartClick Edge to the side of the device

4. Once both SmartClick Edges are mounted, you can attach the SmartClick Rail(s) ⑤. To do this, hold the SmartClick Rail at a 45-degrees angle (with the top facing the device), insert it into the recess at the top, push it upward, then tilt the bottom side toward the device and press the SmartClick Rail downward.

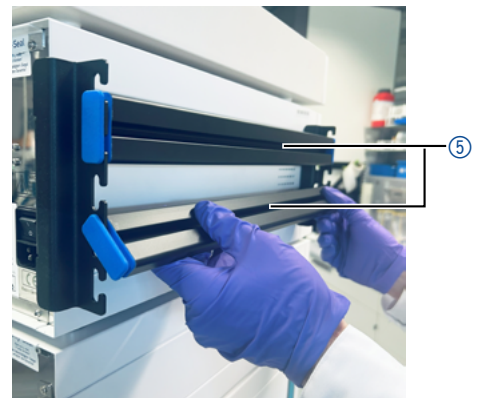


Fig.4: Attaching the SmartClick Edge to the side of the device

Result The SmartClick Edges and the SmartClick Rail(s) have been installed and are now ready for use.

Next steps Install the desired components on the SmartClick Rail(s).



Note: When attaching the desired components, tilt the SmartClick adapter 45 degrees with the top facing the device, insert the adapter into the rail, and press down on the front of the adapter.



Note: Some KNAUER accessories come with AZURA Click pre-installed. First, remove AZURA Click, then install the AZURA SmartClick adapter A700884. Use the screws that were used to install AZURA Click.

For more information about AZURA SmartClick, please visit the KNAUER website:

